## INSTRUCTION MANUAL

### VWR® Water bath W20

**European Catalogue Number:**  
720-2423

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**Gebrauchsanweisung (DE) 10-13**

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Version: 1  
Issued: 22122014

Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

The unit can have hot surfaces and hot water

Do not use the unit in an explosive environment

## Avoiding Electrical Shock

Make sure that the voltage rating on the unit corresponds to the mains power voltage

Make sure the mains power socket (wall socket) has a functioning electrical ground connection 

Make sure the units power plug corresponds to the mains power socket

Always use the units grounded power cord between the unit and mains power socket

Do not use the unit if the power cord is damaged or if the unit is damage

Do not open the unit, all repairs should be done by an authorized technician

---

## Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorisations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorisation, unless any refusal is due to a defect of the product.

## Equipment disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Description	Quantity
Water Bath W20	1
Power cord (EU-plug, UK-plug)	2
Instruction manual	1

## Installation

Unpack the delivery. Check that the unit is not damaged and that all parts are included.

Place the unit on the work area. Connect the unit to an electrical grounded mains power socket, use the units power cord (which has a with ground connection).

Check that the unit is OFF; no lights on the power switch or display. (If the unit is ON, green light on the power switch and red light on the display, then press the down side of the power switch, the lights goes OFF.)

## Intended use and overview

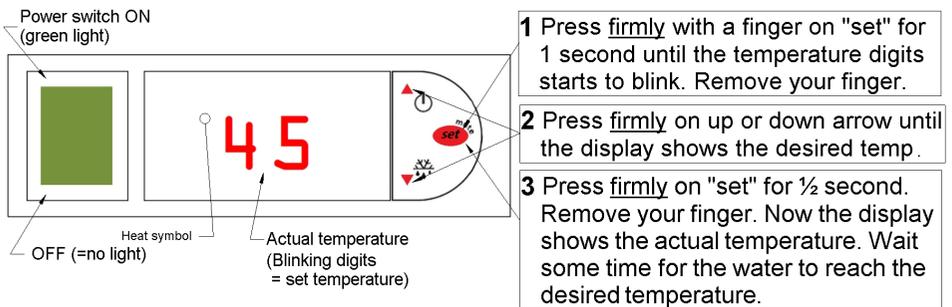
The product is for use at a pathology/histology laboratory for in vitro diagnostics, for preparation of human tissue sections on glass slides for microscopy. The units casing is made of white powder coated aluminium. The water bath (inside) is coated with black corrosion resistant thermoplastic. The unit is equipped with a power switch with green light, a temperature control unit with digital display, internal heating element and an overheat protection fuse.

The flat black surface is heated and can be used for drying and heating samples. The temperature is slightly lower than the water temperature.

An optional lid is available to cover the water. It prevents evaporation and dust contamination when the unit is not in use. VWR order number: 720-2424

## Product Specifications

Mains power:	230 VAC 50/60Hz
Power consumption:	190 W
Temperature range:	+20 to +75°C
Thermostat:	Digital
Temp. accuracy:	±1°C
Water capacity:	1,5 L
External dimensions:	W:247 x D:280 mm x H:78 mm
Internal dimension:	W:198 x D:156 x H:60 mm
Weight:	2,4 kg

## Getting Started

Fill the water bath with 1,3 litres of water (distilled water to avoid deposits).

Turn the unit ON by pressing the upper part of the power switch; the power switch shows green light and the display shows actual water temperature.

To change the desired temperature, see instructions **1-2-3** in the figure above.

If you press longer than 3 seconds and the display shows **P5** then reset the unit by pressing the power switch OFF and then ON. Start at 1 again, see above.

## Important Guidelines

Be careful when the emptying the water bath. First turn the unit OFF. Then disconnect the power cord first from the mains supply and then from the unit. Please observe that the water can be hot. Carefully empty the water from the side of the unit. Do not let water enter the back or front of the unit. Wipe off the unit from any remaining water, leaving a dry surface.

Service and repair should only be made by an authorized technician.

## Accessories

Description	Cat. No.
Lid. Covers the water, prevents evaporation and dust contamination	720-2424

Review the information in the table below to troubleshoot operating problems.

Problem	Cause	Solution
No heat	Too low set temperature	Increase the desired (set) temperature
	Heating element problem	Repair by an authorized technician
No power, no light	Power cord not firmly attached	Check the connection to mains power and to the unit
	The units fuse has blown	Replace the units fuse, located under the power cord socket
	The internal thermo fuse has blown	Repair by an authorized technician
Power switch has green light but no light on the display	Control unit not functioning	Repair by an authorized technician

## Technical service

### Web Resources

Visit the VWR website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to the VWR Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit [www.vwr.com](http://www.vwr.com).

## Warranty

**VWR International** warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

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# Declaration of Conformity

Undersigned representative for the following Manufacturer:

Manufacturer: KUNZ INSTRUMENTS AB  
Address: Tradgardsgatan 9, SE-14934 Nynashamn, Sweden

Declare hereby that the following product: 720-2423 Water bath W20

Comply with the following EC Directive and standards:

Directive	Title
98/79/EC	In Vitro Diagnostic Medical Devices (IVD), Annex III: other
2011/65/EU	RoHS2

Standard	Title
IEC/SS-EN 61010-1:2001	Safety requirements for electrical equipment for measurement, control and laboratory use Part 1: General requirements
IEC/SS-EN 61010-2-101:2002	Safety requirements for electrical equipment for measurement, control and laboratory use Part 2-101: Particular requirements for in vitro diagnostic (IVD) medical equipment
IEC/SS-EN 61326-1:1997	Electrical equipment for measurement, control and laboratory use - EMC requirements
IEC/SS-EN 61326-2-6:2006	Electrical equipment for measurement, control and laboratory use - EMC requirements Part 2-6: Particular requirements - In-vitro diagnostic (IVD) medical equipment
SS-EN ISO 14971:2012	Medical devices - Application of risk management to medical devices

Nynashamn December 4, 2014

---

Anders Olsson, Manufacturers representative

# Local VWR offices in Europe and Asia Pacific

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## India

VWR Lab Products Private Limited  
135/12, Brigade Towers, 2nd Floor  
Front wing, Brigade Road,  
Bengaluru, India – 560 025  
Tel: +91-2522-647911/922 (Mumbai)  
Tel: +91-80-41117125/26 (Bangalore)  
Fax +91-80-41117120  
E-mail: vwr\_india@vwr.com

## Ireland / Northern Ireland

VWR International Ltd / VWR International  
(Northern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: 01 88 22 222  
Fax: 01 88 22 333  
E-mail sales@ie.vwr.com

## Italy

VWR International PBI S.r.l.  
Via San Giusto 85  
20153 Milano (MI)  
Tel.: 02-3320311/02-487791  
Fax: 02-332031307/02-40090010  
E-mail: info@it.vwr.com

## The Netherlands

VWR International B.V.  
Postbus 8198  
1005 AD Amsterdam  
Tel.: 020 4808 400  
Fax: 020 4808 480  
E-mail: info@nl.vwr.com

## New Zealand

VWR International LP  
241 Bush Road  
Albany  
0632 New Zealand

## Norway

VWR International AS  
Haavard Martinsens vei 30  
0978 Oslo  
Tel.: 02290  
Fax: 815 00 940  
E-mail: info@no.vwr.com

## Poland

VWR International Sp. z o.o.  
Limbowa 5  
80-175 Gdansk  
Tel.: 058 32 38 210  
Fax: 058 32 38 205  
E-mail: labart@pl.vwr.com

## Portugal

VWR International - Material de  
Laboratório, Lda  
Edifício Neopark  
Av. Tomás Ribeiro, 43-3 D  
2790-221 Carnaxide  
Tel.: 21 3600 770  
Fax: 21 3600 798/9  
E-mail: info@pt.vwr.com

## Singapore

VWR Singapore Pte Ltd  
18 Gul Drive  
Singapore 629468  
Tel: +65 6505 0760  
Fax: +65 6264 3780  
E-mail: sales@sg.vwr.com

## Spain

VWR International Eurolab S.L.  
C/ Tecnología 5-17  
A-7 Llinars Park  
08450 - Llinars del Vallès  
Barcelona  
Tel.: 902 222 897  
Fax: 902 430 657  
E-mail: info@es.vwr.com

## Sweden

VWR International AB  
Fagerstagatan 18a  
163 94 Stockholm  
Tel.: 08 621 34 00  
Fax: 08 621 34 66  
E-mail: info@se.vwr.com

## Switzerland

VWR International GmbH  
Lerzenstrasse 16/18  
8953 Dietikon  
Tel.: 044 745 13 13  
Fax: 044 745 13 10  
E-mail: info@ch.vwr.com

## Turkey

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.  
a VWR International Company  
Orta Mah. Cemal Gürsel Caddesi  
Ördekcioglu İşmerkezi No.32/1  
34896 Pendik - İstanbul  
Tel.: +90216 598 2900  
Fax: +90216 598 2907  
Email: info@pro-lab.com.tr

## UK

VWR International Ltd  
Customer Service Centre  
Hunter Boulevard  
Magna Park  
Lutterworth  
Leicestershire  
LE17 4XN  
Tel.: 0800 22 33 44  
Fax: 01455 55 85 86  
E-mail: uksales@uk.vwr.com