

**SOFTWARE SETUP
INSTALLAZIONE SOFTWARE
INSTALACIÓN SOFTWARE
CONFIGURATION DU LOGICIEL
SOFTWARE EINSTELLUNG
INSTALAÇÃO SOFTWARE**

Model
BL224T1
BL254T1

Version: 1
Issued: 17, 12, 2015



1. Start the Software

Double click on the icon.

1. Aprire il Software

Cliccare due volte sull'icona.

1. Iniciar Software

Doble click sobre el icono.

1. Lancez le logiciel


Cliquez deux fois sur l'icône.

1. Software Starten

Doppelklick auf die Ikone.

1. Iniciar Software

Doble click sobre o ícone.

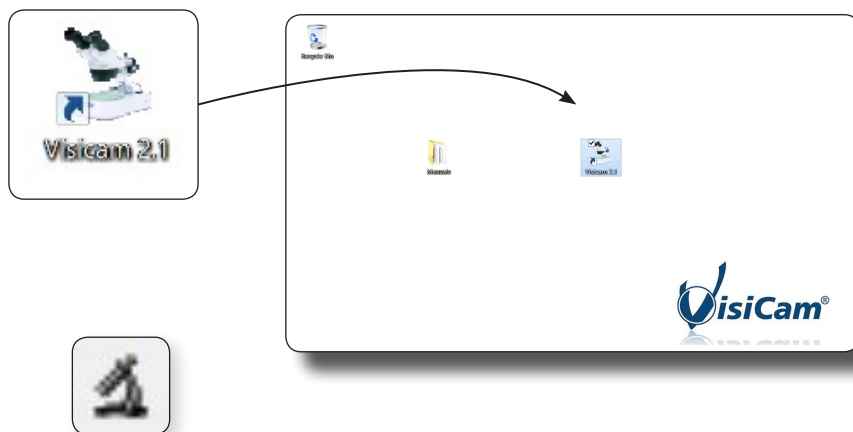

2. How to set the live image

Press the icon.

2. Come impostare le immagini in live

Cliccare sull'icona.

2. Como ver imagen en tiempo real

Click en el icono.

2. Comment régler l'image en direct


Appuyez sur l'icône.

2. Einstellung der live image


Drücken Sie die Ikone.

2. Como ver imagem em tempo real

Click em o ícone.

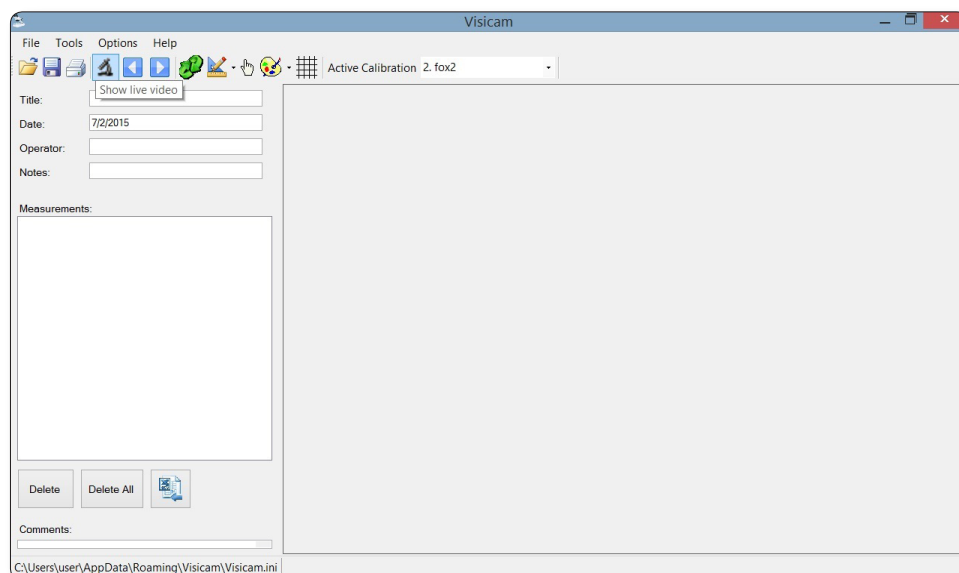

Select the device "Camera-X3"
Selezionare il dispositivo "Camera-X3"
Seleccionar la unidad "Camera-X3"
Sélectionnez l'appareil "Camera-X3"
Wählen Sie "Camera-X3" aus
Selecionar a unidade "Camera-X3"

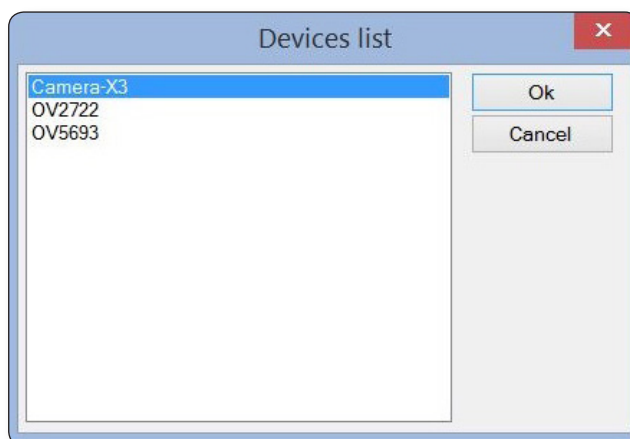

3. How to set the colors parameter

Set the screen brightness at the maximum.

3. Come impostare i parametri di colore

Impostare al massimo la luminosità dello schermo.

3. Como configurar los parámetros de color

Posicionar el brillo de la pantalla al máximo.

3. Comment définir le paramètre de couleurs


Régler la luminosité de l'écran au maximum.

3. Einstellung der Farbparameter

Stellen Sie den Bildschirm auf die maximale Helligkeitseinstellung ein.

3. Como configurar os parâmetros de cor

Posicionar o brilho de la tela al máximo.

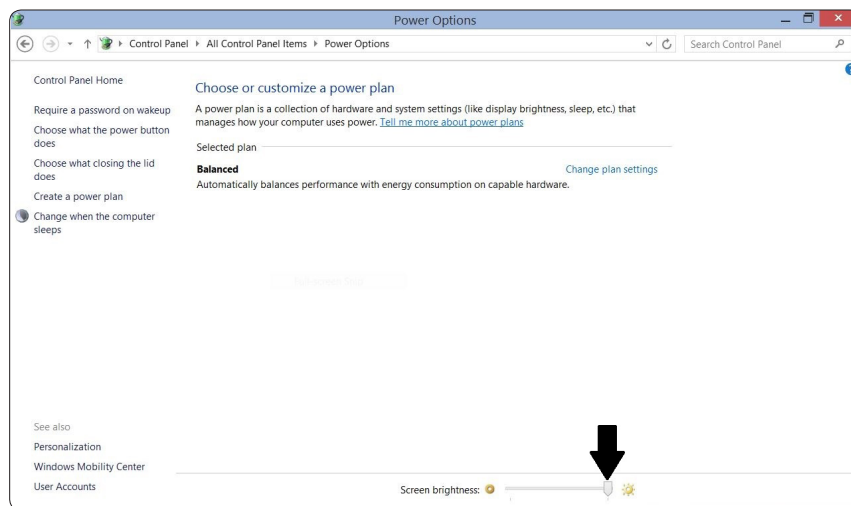

Open the settings of the software (icon)

Set the contrast value from 9 to 12 as needed.

Aprire il menu delle impostazioni del software (icona).

Impotare il valore del contrasto tra 9 e 12 in base alla necessità.

Ouvrez les paramètres du logiciel (icône)

Réglez la valeur du contraste du 9 au 12 comme requis.

Abrir configuración de software (icono)


Posicionar el valor de contraste entre 9 y 12 según sus necesidades.

Öffnen Sie die Software Einstellungen (Ikon)

Stellen Sie den Kontrastwert von 9 bis 12 ein, wie benötigt.

Abrir configuração do software (Ikon)

Posicionar o valor do contraste entre 9 y 12 segun as necessidades.

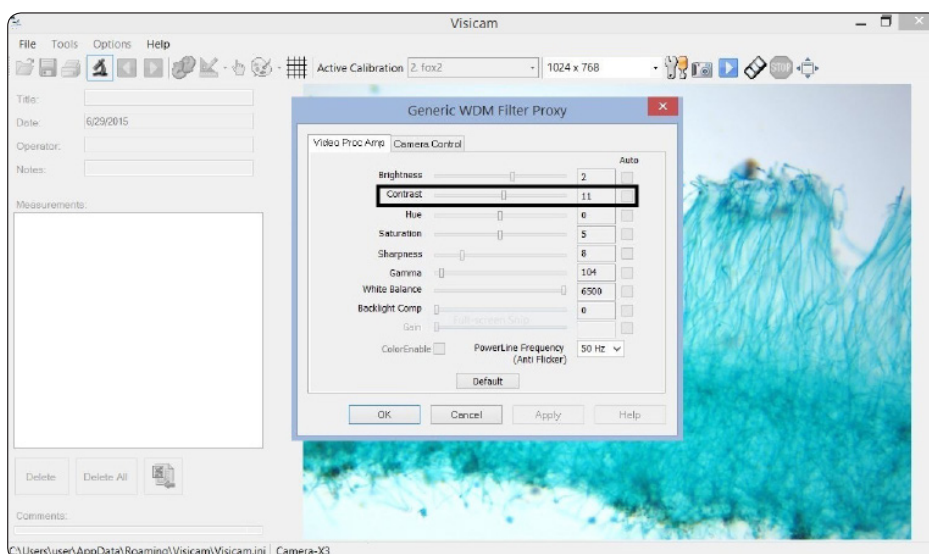

Alemanha

VWR International GmbH
Hilpertstraße 20a
D - 64295 Darmstadt
Freecall: 0800 702 00 07
Fax: 0180 570 22 22*
Email: info@de.vwr.com
*0,14 €/Min. aus d. dt. Festnetz

Austrália

VWR International, Pty Ltd.
1/31 Archimedes Place
Murarrie, Queensland, 4172
Tel.: 1300 727 696
Fax: 1300 135 123

Áustria

VWR International GmbH
Graumannsgasse 7
1150 Wien
Tel.: 01 97 002 0
Fax: 01 97 002 600
E-mail: info@at.vwr.com

Bélgica

VWR International bvba
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: 016 385 011
Fax: 016 385 385
E-mail: customerservice@be.vwr.com

China

VWR International China Co., Ltd
Rm.219, 2100 Dongming Road
Pudong New District
Shanghai 200123
Tel.: +86-21-5898 6888
Fax: +86-21-5855 8801
E-mail: info_china@vwr.com

Dinamarca

VWR - Bie & Berntsen
Transformervej 8
2730 Herlev
Tel.: 43 86 87 88
Fax: 43 86 87 90
E-mail: info@dk.vwr.com

Espanha

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Finlândia

VWR International Oy
Valimotie 9
00380 Helsinki
Tel.: 09 80 45 51
Fax: 09 80 45 52 00
E-mail: info@fi.vwr.com

França

VWR International S.A.S.
Le Périgares – Bâtiment B
201, rue Carnot
94126 Fontenay-sous-Bois cedex
Tel.: 0 825 02 30 30 (0,15 € TTC/min)
Fax: 0 825 02 30 35 (0,15 € TTC/min)
E-mail: info@fr.vwr.com

Índia

VWR Lab Products Private Limited
135/12, Brigade Towers, 2nd Floor
Front wing, Brigade Road,
Bengaluru, India – 560 025
Tel.: +91-80-41117125/26 (Bengaluru)
Tel.: +91-2522-647911/922 (Mumbai)
Fax: +91-80-41117120
E-mail: vwr_india@vwr.com

Hungria

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: (52) 521-130
Fax: (52) 470-069
E-mail: info@hu.vwr.com

Irlanda / Irlanda do Norte

VWR International Ltd /
VWR International (Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Itália

VWR International PBI S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 800-152999/02-40090010
E-mail: info@it.vwr.com

Noruega

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Nova Zelândia

Global Science - A VWR Company
241 Bush Road
Albany 0632, Auckland
Tel.: 0800 734 100
Fax: 0800 999 002
E-mail: sales@globalscience.co.nz

Países Baixos

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

Polónia

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: 58 32 38 200 do 204
Fax: 58 32 38 205
E-mail: info@pl.vwr.com

Portugal

VWR International - Material de Laboratório, Lda
Edifício Neopark
Av. Tomás Ribeiro, 43- 3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Reino Unido

VWR International Ltd
Customer Service Centre
Hunter Boulevard - Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

República Checa

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Stříbrná Skalice
Tel.: +420 321 570 321
Fax: +420 321 570 320
E-mail: info@cz.vwr.com

Singapura

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Suécia

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: kundservice@se.vwr.com

Suiça

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

Turquia

Pro Lab Laboratuvar Teknolojileri Ltd.Şti.
a VWR International Company
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - İstanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info@pro-lab.com.tr

**VISITE-NOS EM WWW.VWR.COM
PARA CONHECER AS ÚLTIMAS
NOVIDADES, OFERTAS ESPECIAIS
E INFORMAÇÕES DETALHADAS**
