



HistoChoice® Mounting Media

Code	Description	Size
H157-120ML	HistoChoice® Mounting Media	120 mL
H157-475ML	HistoChoice® Mounting Media	475 mL

General Information

HistoChoice® Mounting Media was designed for use with HistoChoice® Clearing Agent, but can also be used as a separate, stand-alone medium for standard slide preparation. It is suitable for both thick and thin specimen mounting and is compatible with immersion oil. The low viscosity of this mounting medium facilitates easy application and bubble-free preparations, which saves valuable lab time.

- Does not crack or splinter with age and will not discolor
- Low viscosity for bubble-free slide preparations
- Works for all types/sizes of specimens

Storage/Stability

Store at room temperature (18 to 26°C).

Product Use Limitations

For research use only. Not for therapeutic or diagnostic use.

Protocol/Procedure

Mounting specimens with HistoChoice® Mounting Media

1. Drain excess clearing agent from the slide that will be coverslipped.
2. Lay the slide on a level surface and gently apply a drop or two of HistoChoice® Mounting Media to cover the specimen.
3. Apply a coverslip to the slide by lowering it slowly from a 45° angle until it covers the specimen and the mounting media has spread under the glass.
4. Allow sample to dry.

Frequently Asked Questions

What is the best way to remove a coverslip from a specimen?

- To remove a coverslip mounted with HistoChoice® Mounting Media, soak the slide in HistoChoice® Clearing Agent for 1 – 2 hours or dip in a Coplin jar of xylene.

For Technical Support

Toll Free: 1-800-610-2789 (USA & Canada)

Fax: (440) 349-0235

Email: techinquiry@amresco-inc.com

AMRESKO, LLC

A VWR Company

Corporate Headquarters
28600 Fountain Parkway
Solon, Ohio USA 44139-4300

Tel: 440/349-1199

Fax: 440/349-1182

www.amresco-inc.com

HistoChoice® Mounting Media

ZY0588

Rev. 1 12/2015

© Copyright 2010 by AMRESKO, LLC

All Rights Reserved.

AMRESKO® is a registered trademark of AMRESKO, LLC