EN - English		1
FR - Français		13
ES - Español		26
IT - Italiano		39
DE - Deutsch		52
PT - Português		65
NL - Nederlands		69
NO - Norsk		73
DA - Dansk		77
SV - Svenska		81
FI - Suomi		85
HU - Magyar		89
PL - Polski		93

Instruction Manual

Digital & Analogue

Heavy Duty Vortex Mixers

Microplate Vortex Mixers

European Catalogue Numbers:

Analogue Heavy Duty Vortex Mixers

Euro plug and Swiss adapter: 444-7133

UK plug: 444-7134

Digital Heavy Duty Vortex Mixers

Euro plug and Swiss adapter: 444-7130

UK plug: 444-7131

Analogue Microplate Vortex Mixers

Euro plug and Swiss adapter: 444-0986

UK plug: 444-0993

Digital Microplate Vortex Mixers

Euro plug and Swiss adapter: 444-0985

UK plug: 444-0992

TABLE OF CONTENTS

Package Contents	1
Warranty	1
Installation	2
Maintenance & Servicing	2
Environmental Conditions	2
Safety Instructions	3
Warning	3
Standards & Regulations	3
Specifications	4
Vortex Mixer Control Panels	4-5
Introduction	6
Operating Instructions	6-7
Head/Accessory Installation Instructions	8-9
Troubleshooting	10
Accessories	11-12
EC Declaration of Conformity	97-98

PACKAGE CONTENTS

Heavy Duty Vortex Mixer,	OR	Microplate Vortex Mixer,
Analogue or Digital		Analogue or Digital
183cm detachable power cord		183cm detachable power cord
with Euro plug and UK plug or Swiss adapter*		with Euro plug and UK plug or
(1) cup head		Swiss adapter*
(1) universal holder		(1) cup head
(1) universal holder cover		(1) microplate holder, single
(1) foam insert (1,5-2,0 mL microtubes, holds 38)		Instruction manual
Instruction manual		

**Dependent upon model*

WARRANTY

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of delivery. If a defect is present, VWR will, at its option and cost, repair, replace, or refund the purchase price of this product to the customer, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. If the required maintenance and inspection services are not performed according to the manuals and any local regulations, such warranty turns invalid, except to the extent, the defect of the product is not due to such non-performance.

Items being returned must be insured by the customer against possible damage or loss. This warranty shall be limited to the aforementioned remedies. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Compliance with local laws and regulations

The customer is responsible for applying for and obtaining the necessary regulatory approvals or other authorizations necessary to run or use the Product in its local environment. VWR will not be held liable for any related omission or for not obtaining the required approval or authorization, unless any refusal is due to a defect of the product.

INSTALLATION

Upon receiving the VWR Vortex Mixer, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, remove the protective coverings from the feet and place the Vortex Mixer on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is clean and free of dust. Always place the unit on a sturdy work surface.

The Vortex Mixer is supplied with a power cord that should be plugged into a properly grounded outlet. If the supplied power cord does not meet your needs, please use an approved power cord that suits local codes and electric supply. Replacement of the plug must be made by a qualified electrician.

MAINTENANCE & SERVICING

The Vortex Mixer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly. **DO NOT** use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

ENVIRONMENTAL CONDITIONS

Operating Conditions: Indoor use only.

Temperature: 18 to 33°C

Humidity: 20% to 85% relative humidity, non-condensing

Altitude: 0 to 2000 M above sea level

Non-Operating Storage:

Temperature: -20 to 65°C

Humidity: 20% to 85% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

EQUIPMENT DISPOSAL

This equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment at life-cycle end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical, and/or radiological contamination, so as to protect the persons involved in the disposal and recycling of the equipment from health hazards. For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

SAFETY INSTRUCTIONS

Please read the entire instruction manual before operating the Vortex Mixer.

WARNING! DO NOT use the Vortex Mixer in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift the Vortex Mixer by the head. All heads, including cup head, are removable. They can pop off if lifted by them.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

STANDARDS & REGULATIONS

VWR international hereby declares under its sole responsibility that the products comply with the requirements of the following directives and associated standards:

Associated EU guidelines:

EMC directive 2014/30/EU

LVD directive 2014/35/EU

RoHS directive 2011/65/EU

Safety standards:

IEC 61010-1

Safety requirements for electrical equipment for measurement, control and laboratory use. Part I: General Requirements.

IEC 61010-2-051

Part II: Particular requirements for laboratory equipment for mixing and stirring.

UL Std. No. 61010-1

EMC standards:

EN61326-1

FCC Part 15 Subpart B

HEAVY DUTY VORTEX MIXER SPECIFICATIONS

Overall dimensions (L x W x H):	24,1 x 16,8 x 16,5cm
Electrical (50/60 Hz):	0,13 amps 30 watts
Fuses:	5mm x 20mm, 5 amp 250 volt quick acting
Speed range	On Mode: 300-2500rpm
	Touch Mode: 300-3500rpm
Orbit:	4,9mm
Duty rating:	Continuous duty
Controls:	Analogue: Touch/off/on - rocker switch
	Speed knob: variable 1 to 10 dial marks
	Digital: See page 5
Capacity:	1,1kg
Ship weight:	6,8kg
European Catalogue Numbers:	
	Analogue: Euro plug and Swiss adapter: 444-7133
	UK plug: 444-7134
	Digital: Euro plug and Swiss adapter: 444-7130
	UK plug: 444-7131

MICROPLATE VORTEX MIXER SPECIFICATIONS

Overall dimensions (L x W x H):	26,6 x 13,7 x 11,4cm
Electrical (50/60 Hz):	0,13 amps 30 watts
Fuses:	5mm x 20mm, 5 amp 250 volt quick acting
Speed range	On Mode: 300-2500rpm
	Touch Mode: 300-3500rpm
Orbit:	3,5mm
Duty rating:	Continuous duty
Controls:	Analogue: Touch/off/on - rocker switch
	Speed knob: variable 1 to 10 dial marks
	Digital: See page 5
Capacity:	2 Microplates
Ship weight:	5,4kg
European Catalogue Numbers:	
	Analogue: Euro plug and Swiss adapter: 444-0986
	UK plug: 444-0993
	Digital: Euro plug and Swiss adapter: 444-0985
	UK plug: 444-0992

ANALOGUE VORTEX MIXER CONTROL PANEL

The front panel of the Analogue Vortex Mixer contains all the switches and control knob needed to operate the unit.

A. 3-way rocker switch: touch/off/on rocker switch starts/stops the vortexing function.

B. Speed control: The speed control knob controls the vortexing speed. User should turn knob clockwise until the vortexer reaches the desired speed.

DIGITAL VORTEX MIXER CONTROL PANEL

The front panel of the Digital Vortex Mixer contains all the switches, controls and displays needed to operate the unit.

A. 3-way rocker switch: touch/standby/on rocker switch starts/stops the vortexing function.

B. Time display: Displays accumulated time (continuous mode) or how much time is remaining (timed mode). The display range is from 0 to 9,999 minutes in one (1) second increments. The display will indicate minutes and seconds until the timer reaches 99 minutes and 59 seconds (99:59), then the display will automatically display minutes up to 9,999.

C. Up/down arrows for set-point control.

D. Speed display: Displays the speed of the Vortex Mixer.

E. Up/down arrows for set-point control.

VORTEX MIXER INTRODUCTION

Your Vortex Mixer is ready for most one-handed applications. The Vortex Mixer operates by mixing samples just prior to testing.

TOUCH MODE OPERATION

NOTE: Due to the robust mechanical design of this unit, more downward force is required to operate this unit in Touch Mode to mix single tubes as compared to other models. Another option is to set your unit to the ON mode for mixing a single tube.

TOUCH MODE OPERATION

To ensure the stability of your vortex mixer, it is very important to start with a smooth, clean work surface and the mixers (4) feet. If the mixer starts to “walk”, remove the power to the unit and clean the work surface and the (4) feet with isopropyl alcohol. Press the unit down firmly on the work surface to ensure the feet are properly bonded.

ANALOGUE VORTEX MIXER OPERATING INSTRUCTIONS

1. Make sure the 3-way rocker switch is in the center, off position. Plug the power cord into a properly grounded outlet.
2. For continuous operation, push the rocker switch to the right, on position. For intermittent/touch operation, push the rocker switch to the left, touch position.
3. In either case, turn the speed knob to the desired setting. In the on position, you will see the head in motion immediately. In the touch position, the vessel you are using must be pushed down on the head attachment to achieve motion. To get the desired mix, vary the speed, using the speed knob, and/or vary the angle of contact and pressure against the head.
4. When finished with either continuous or intermittent/touch operation, return rocker switch to the center, off position.
5. The continuous, on mode is for the accessories.
6. The intermittent, touch mode is intended for short mixing times of one (1) minute or less at full speed.
7. To completely cut off power to the unit, disconnect the power cord from the wall outlet.

DIGITAL VORTEX MIXER OPERATING INSTRUCTIONS

1. Make sure the 3-way rocker switch is in the center, standby position. Plug the power cord into a properly grounded outlet. The unit is now on and in a standby mode.
2. For continuous operation, push the rocker switch to the right, on position. For intermittent/touch operation, push the rocker switch to the left, touch position.
3. In either case, set the speed. In the on position, you will see the head in motion immediately. In the touch position, the vessel you are using must be pushed down on the head attachment to achieve motion. To get the desired mix, vary the speed, using the speed up/down arrows, and/or vary the angle of contact and pressure against the head.

Setting speed:

To set the speed press the up/down arrows below the speed display until you reach the desired speed.

- Speed while running in the on position is 300 to 2500rpm
- Speed while running in the touch position is 300 to 3500rpm

Setting time:

Time operation works in both touch and on positions.

- To run in timed mode (programmed time), start by pressing the rocker switch to the left, touch position. Press the up/down arrows below the time display until you reach the desired time. While running unit in timed mode, the time display will show time remaining, counting down to zero (0:00).
- To run in continuous mode (accumulated time), reset the timer to zero (0:00) before running. Time display will show the accumulated time and run until you press the rocker switch to the center, standby position.

4. When finished with either continuous or intermittent/touch mode, return rocker switch to the center, standby position.
5. The continuous, on mode is for accessories.
6. The intermittent, touch mode is intended for short mixing times of one (1) minute or less at full speed.
7. To completely cut off power to the unit, disconnect the power cord from the wall outlet.

Beeper Reference:

To silence the beeper operation (except for error codes), with the unit in standby mode, remove accessories or cup head. Press and hold the head mount. Press the time down arrow. Release the time down arrow. Release the head mount. To restore to factory setting, repeat procedure.

ACCESSORY INSTALLATION INSTRUCTIONS (ALL UNITS)

Installing the cup head, microplate trays, universal holder or other head attachments:

1. Move the 3-way rocker switch to the center, off/standby position.
 2. Place accessory on head mount of the unit, aligning the square on the bottom of the accessory with the square on the top of the unit. Lower accessory until it is sitting on the head mount.
 3. Turn the accessory counter-clockwise until it snaps into place (approximately 1/8 of a turn). Add samples to attachments after the accessory is secured onto the Vortex Mixer. Be sure samples are secure before turning unit on.
 4. To remove an installed accessory, move the 3-way rocker switch to the center, off/standby position. Turn the accessory clockwise (approximately 1/8 of a turn) and then lift the accessory straight up. Remove samples before removing the accessories.
- To load a microplate into the microplate tray, position the microplate against the foam bumpers first, then push down the other side to ensure a secure fit. To release a microplate, push it against the microplate tray's foam bumper and lift to remove.
 - Maximum speed may vary when vortexing microplates. Speed may be dependant on the height of the skirt on the microplate or material of the microplate.
 - Rubber vibration dampeners on the bottom of the microplate tray and universal holder may need to be reinserted into the proper position when vortexing at high speeds. Squeeze firmly as you maneuver it back into position.

ACCESSORY INSTALLATION INSTRUCTIONS (HEAVY DUTY VORTEX MIXER)

Installing the vessel holders:

1. Attach the Universal Holder to the vortex mixer.
2. Place the Grip Mat in the bottom of the Universal Holder.
3. Place the flask on top of the Grip Mat, in the center of the holder.
4. Pull the center hole of the Vessel Holder over the sample container.
5. With the arrow on the Vessel Holder close to the front of the unit, the straps should line up with the posts of the Universal Holder.
6. Push the hole on the strap over the posts on the holder. Secure the strap so that it fits into the slot on the holder. Repeat this for the remaining two straps.
7. Be sure the sample is secure before turning unit on.


Installing the Universal Holder Cover:

1. Place the Foam Insert into the Universal Holder.
2. Lay the Universal Holder Cover on top of the foam, and align the edges.
3. Push the hole on the cover down over the post on the holder. Squeeze the strap so that it fits into the slot on the holder, leaving the cover sitting flat on the foam. Repeat this process for the remaining two posts.
4. To remove the accessory cover, secure the straps and pull straight up.

ACCESSORY INSTALLATION INSTRUCTIONS (HEAVY DUTY VORTEX MIXER)

Installing Single Tube Holder:

1. Move the 3-way rocker switch to the center, off/standby position.
2. Remove cup head, or any other accessory from the Vortex Mixer. Turn the cup head clockwise, approximately 1/8 of a turn, (see figure 1) and then lift the accessory straight up (see figure 2). Remove samples before removing the accessories.
3. Align the base plate of the Single Tube Holder over the head mount of the Vortex Mixer with the two Single Tube Holder alignment posts at the back of the unit. Lower Single Tube Holder onto Vortex Mixer, (see figure 3) the magnets will hold the Single Tube Holder in place. To realign this attachment, pull up from the handle until the magnets release, check the alignment and place again.
4. Now that the Single Tube Holder is in place, attach the cup head on the head mount of the unit. Turn the cup head counter-clockwise until it snaps into place, approximately 1/8 of a turn. (see figure 4)
5. Lift up on the spring loaded upper portion of the Single Tube Holder and slide a 6,4 to 11,4cm long tube on the cup head so that the tube is aligned in the middle (see figure 5). You are now ready to mix your sample.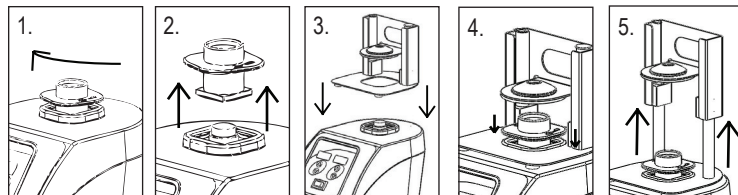

TECHNICAL SERVICE

For more information or technical service contact your local VWR representative or visit www.vwr.com.


Web resources: Visit VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional products information and special offers

TROUBLESHOOTING (ALL UNITS)

Error	Cause of Error	How to Fix
Analogue Models Motor does not run in the touch or on mode Digital Models E3 Error	Analogue and Digital Models There is no motion on the motor or the motor is not working properly	Analogue and Digital Models 1. Turn the rocker switch to the center off/standby position 2. Check to verify the load has not been exceeded 3. Check and clear any possible obstructions of the head mount 4. If the unit still does not work, please contact your VWR representative

ACCESSORIES FOR THE HEAVY DUTY VORTEX MIXER

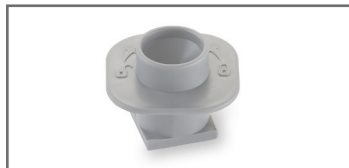Cup Head

Replacement for cup head supplied with Heavy Duty Vortex Mixer. Designed for mixing one tube at a time.

Cat. No. 444-7136

Universal Holder

Replacement for Universal Holder supplied with Heavy Duty Vortex Mixer to mix irregularly shaped objects when Universal Holder Cover is in place. Also used as a holder for the foam tube inserts. Measures 15,5x15,0x7,6cm.

Cat. No. 444-7139

Universal Holder Cover

Replacement for Universal Holder Cover supplied with Heavy Duty Vortex Mixer. The cover allows for mixing irregularly shaped objects.

Cat. No. 444-7140

Single Tube Holder

Single tube hands free mixing designed to fit on the VWR Heavy Duty Vortex Mixer. Easily attached to the top of mixer and is magnetically secured. Accepts tubes from 6,4-11,4cm tall. Minimum tube diameter is 19mm.

Cat. No. 444-7150

Microplate Holder (single)

Designed to hold one microplate.

Cat. No. 444-7137

Microplate Holder (double)

Designed to hold two standard microplates.


Cat. No. 444-7138

Microplate Holder (quad)

Designed to hold four standard microplates.

Cat. No. 444-7151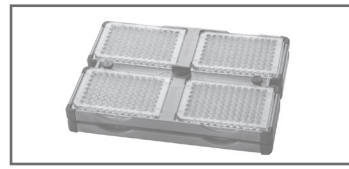

Stackable Microplate Holder Four

Designed to maximize the capacity of the Heavy Duty Vortex Mixer to eight microplates by stacking the tray on top of the four plate holder.

Cat. No. 444-7152

ACCESSORIES FOR THE HEAVY DUTY VORTEX MIXER

Foam Insert for 0,5mL microtubes

Foam insert holds (52) 0,5mL microtubes.

Cat. No.

444-7141

Foam Insert for 1,5 to 2,0mL microtubes

Foam insert holds (38) 1,5 to 2,0mL microtubes.

Cat. No.

444-7142

Foam Insert for 12-13mm test tubes

Foam insert holds (34) 12-13mm diameter test tubes.

Cat. No.

444-7143

Foam Insert for 15-18mm test tubes

Foam insert holds (20) 15-18mm diameter test tubes. Ideal for 15mL centrifuge tubes.

Cat. No.

444-7144


Foam Insert for 19-21mm test tubes

Foam insert holds (18) 19-21mm diameter test tubes.

Cat. No.

444-7145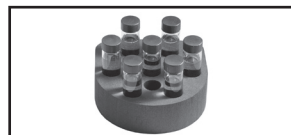


Foam Insert for 22-25mm test tubes

Foam insert holds (13) 22-25mm diameter test tubes.

Cat. No.

444-7146

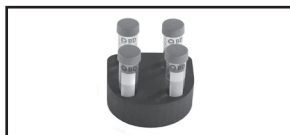Foam Insert for 26-29mm test tubes

Foam insert holds (4) 26-29mm diameter test tubes. Ideal for 50mL centrifuge tubes.

Cat. No.

444-7147

Foam Insert Blank

Foam insert without tube openings.

Cat. No.

444-7153

Small Vessel Holder

Rubber holder secures 125 and 250mL Erlenmeyer flasks. Vessel holder also includes a grip mat.

Cat. No.

444-7148


Large Vessel Holder

Rubber holder secures 500 and 1000mL Erlenmeyer flasks. Vessel holder also includes a grip mat.

Cat. No.

444-7149

ACCESSORIES FOR THE MICROPLATE VORTEX MIXER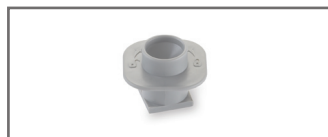


Cup Head

Replacement for cup head supplied with Microplate Vortex Mixer. Designed for mixing one tube at a time.

Cat. No.

444-7136

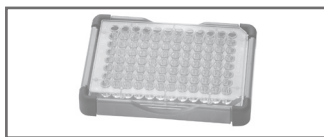Microplate Holder (single)

Replacement for Microplate Holder supplied with Microplate Vortex Mixer. Designed to hold one microplate.

Cat. No.

444-7137

Microplate Holder (double)

Designed to hold two standard microplates.

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The undersigned states that this declaration is issued under the sole responsibility of the manufacturer, VWR International bvba, and that the object(s) of the declaration described in annex 1 are in conformity with the relevant Union harmonization legislation.
In the event of unauthorized modification of any of the products listed in the annex 1, this declaration becomes invalid.

Object Name: VWR Vortex Mixer

Relevant EC Directives: Low Voltage Directive 2014/35/EU
EMC Directive 2014/30/EU
RoHS 2011/65/EU

Harmonized Standards: IEC/EN 61010-1
IEC/EN 61010-2-051
IEC/EN 61326-1
FCC Part 15 Subpart B

Document Number: 14871.2016.0001b

Place & date of issue: Leuven, 25/10/2016

Signed for and on behalf of VWR International bvba:

Yves Van Damme
European Regulatory Affairs Manager

Frankie Vangeel
Managing Director

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India

VWR Lab Products Private Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road, Peenya Post,
Bangalore, India - 560058
Tel.: +91-80-28078400
Fax: +91-80-41117120
Email: vwr_india@vwr.com

Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
Email sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: 02-3320311/02-487791
Fax: 02-332031307/02-40090010
Email: info.it@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
Email: info.nl@vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 22 90 00 00
Fax: 815 00 940
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbow 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax: 058 32 38 205
Email: info.pl@vwr.com

Portugal

VWR International - Material de Laboratório, Lda
Centro Empresarial de Alfragide
Rua da Indústria, nº 6
2610-088 Alfragide
Tel.: 21 3600 770
Fax: 21 3600 798/9
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Linars Park
08450 - Linars del Vallès
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
Email: info.ch@vwr.com

Turkey

VWR International Laboratuvar
Teknolojileri Ltd.Şti.
Orta Mah. Cemal Gürsel Caddesi
Ördekcioglu İşmerkezi No.32/1
34896 Pendik - İstanbul
Tel.: +90216 598 2900
Fax: +90216 598 2907
Email: info.tr@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
Email: uksales@vwr.com

Legal Manufacturer:

VWR International bvba • Researchpark Haasrode 2020 • Geldenaaksebaan 464 • 3001 Leuven • + 32 16 385011 • <http://be.vwr.com>
715124-00 (Rev 5)